



# **INSTANT MILK MICROBIOLOGICAL TEST KIT**







## INSTANT MILK MICROBIOLOGICAL TEST KIT


### SALIENT FEATURES:

- Reveal masking of spoiled milk
- Helps farmer in grading of raw milk for further processing
- No special Laboratory is required
- Result in 10 minutes

### TEST PROCEDURE:

- Set incubator at 37 °C
- Take out microwell from plate and put in incubator
- Take 200ul milk sample into microwell
- Add 20ul reagent solution and mix well
- Incubate at 37 °C for 10 minutes
- Compare color of milk sample with color chart and interpret the result

### RESULT INTERPRETATION:

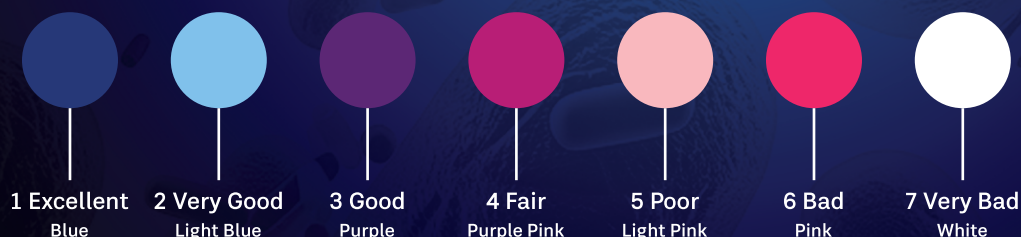

### KIT STORAGE TEMPERATURE:

Store at room temperature

### ORDERING INFORMATION:

CAT. NO.	DESCRIPTION	PACKAGING SIZE
RB2302DD	INSTANT TEST KIT	192 TESTS

### KIT COMPONENTS:

- Microwell plates... 96x2
- Reagent...4ml
- Kit insert

### **Required but not Provided:**

- Micropipette
- Mini incubator

