



Hydrogen Peroxide Rapid Test Strip For Milk & Water

1. Sample:

Raw Milk, Pasteurized Milk, UHT Milk & Water.

2. Operation:

Soak the White Reaction Area of the Strip Into The Sample, After 2 Sec, Take Out.

3. Read The Result:

After 30 Sec, compare the color of the strip with the Following picture (other time is invalid)

Raw	0	4	5	7	8	10
Finished	0	2	5	10	15	20
Water	0	0.5	1	2	5	10

Standard color Scale: mg / L

Notice:

1. Lab Temperature: 15°C - 35°C Relative Humidity \leq 75%
2. The strip is highly susceptible to moisture, light and heat. Please use immediately After take out the strip from the bottle. Cover the rest tightly in bottle.
3. Don't touch the reaction area of the strip. One strip can't be use for twice. The desiccant Can't be take out.

Storage:

2°C - 8°C in cool & Dry Place.
Expiry 12 Months

exons.pk@gmail.com | www.exonscientific.pk

HEAD OFFICE: OFFICE # 215-216, MAMOONA BUILDING, FARID COURT ROAD, LAHORE, PAKISTAN. TEL: +92 42 37 23 2519