



Neutralizers(alkaline Matter) Rapid Test Strip For Milk

1. Sample:

Raw Milk

2. Operation:

Soak the Yellow Reaction Area of the Strip Into The Milk, After 2 Sec, Take Out.

3. Read The Result:

After 1 min, compare the color of the strip with the Following picture (other time is invalid)



Notice:

- 1. Lab Temperature: 15°C - 35°C Relative Humidity \leq 75%**
- 2. The Strip Is Highly Susceptible To Moisture, Light And Heat. Please Use Immediately After Take Out The Strip From The Bottle. Cover The Rest Tightly In Bottle.**
- 3. Don't Touch The Reaction Area Of The Strip. One Strip Can't Be Use For Twice. The Desiccant Can't Be Take Out.**

Storage:

**Room Tamperature & Dry Place.
Expiry 12 Months**

exons.pk@gmail.com | www.exonscientific.pk

HEAD OFFICE: OFFICE # 215-216, MAMOONA BUILDING, FARID COURT ROAD, LAHORE, PAKISTAN. TEL: +92 42 37 23 2519